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MODEL VALIDITY

This manual is valid for all allectric stand-alone medical refrigerators **except** for models of series LR 63, MK 1800, UMKS 65xx and UMKS 14xx (those models are described in a separate user manual).

SETTING UP OF THE APPLIANCE

- The floor of the location site should be even and horizontal. Unevennesses may be compensated by fitting the front feet of the refrigerator.
- The rear side of the appliance may face the wall directly and firm.

CURRENT AND VOLTAGE

Type of current (Alternating Current) and voltage at the location must match indications of the type plate positioned on the inner left wall.

STARTING OPERATION OF THE APPLIANCE

Connect the main plug. Set the temperature control knob to position „5“.

TEMPERATURE SETTING

The built-in temperature controller keeps the set temperature constant inside the cooling compartment.

Position of temp. controller	1	7	0
Temperature condition	highest temperature	lowest temperature	Appliance switched off

When handling stored goods or after storing fresh products temperature may change for a short time. Since fresh products are commonly warmer or warm air enters the cooling compartment. This temperature change will be shown short-time by the temperature display. The temperature of goods already stored will only be changed insignificantly.

ATTENTION: IMPORTANT NOTE!

With temperature setting from controller position „7“ on temperature may drop below 0°C!

DEFROSTING

The appliance defrosts automatically. Thaw water will be collected in an evaporator box and evaporates automatically.

SWITCHING OFF

In order to discontinue operation of the appliance for a longer periode of time disconnect the plug or unscrew/switch off corresponding circuit fuses. Forthmore the door of the appliance must remain opened during non-operating time. This will avoid build-up of odour.

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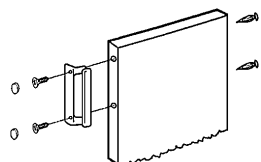
MAINTENANCE AND MALFUNCTION

The appliance doesn't require any maintenance. It was constructed and produced in a way to avoid breakdowns and secure a long lifetime.

Yet in case of troubles occurring during operation please contact us. Quote **in any case the complete model name, spare part index and appliance number**. The type plate giving you this information is positioned on the inner left wall.

You may also establish email contact directly through our homepage www.allectric.com.

ACCESSORIES

The appliance is equipped with light controls including the following functions:

RED LIGHT CONTROL lights up when the inside temperature of the device reaches approximately **+13°C** or more (May also occur with door ajar too long during handling or when temperature controller is set to position „1“).

GREEN LIGHT CONTROL lights up as soon as device is connected to power supply.

CHANGING THE DOOR HINGES

- Unscrew hinge while door closed.
- Lift door off to the bottom.
- Unscrew hinge bolt and screw it into the fixing hole next to it.
- Carefully lift off covers from the left side of the appliance and use them for covering the open fixing holes.
- Put door on its hinges and close it.
- Screw hinges to the left side of the appliance (Don't forget washers).
- If the door shouldn't be in one line to the device casing align the door by adjusting the rectangle holes in the hinge angle (make sure that the screws are connected tightly).
- Unscrew left side door handle.
- Carefully lift off dowel of the right side of the door and use it for sealing the open fixing holes at the left side of the door.
- Screw door handle to the right side of the door.

DOOR HANDLE

The door handle has not been standard mounted with this appliance. By this measure packing volume could be reduced by approximately 40%.

Therefore we may ask you to screw the handle to the appliance on your own.